

DETERMINATION OF ANIONIC SURFACTANTS IN WATER BY FLUORIMETRIC METHOD

INTRODUCTION

Anionic surfactants (A-SAS) represent, by volume, the most important group of surfactants used in cleaning products and the oldest and the largest volume anionic surfactant is soap. In concentrations $0.1 \sim 1.0 \text{ mg/L}$ A-SAS become toxic for aquatic environment.

Lumex Instruments provides sensitive and selective fluorimetric method for the measurement of mass concentration of anionic surfactants in water samples using the FLUORAT-02 analyzer.

The method became the Interstate standard for anionic surfactant in water determination; the benefit of the method is that it doesn't require expensive instrumentation and tedious sample preparation.

MEASUREMENT RANGE

Measurement range	Directives & standards for drinking water	MAC (MPL), mg/L
0.025–1.0 mg/L (0.025–1.0 ppm) (natural, drinking, and waste water)	WHO Guidelines for drinking water quality (2011)	-
	Drinking Water Directive 98/83/EC	-
	US EPA National Secondary Drinking Water Regulations	0.5
	GB 5749-2006 Standards for drinking water quality	0.3
	IS 10500:2012 Drinking water – Specification	0.2
	Resolução CONAMA No 396/2008	_
	Código Alimentario Argentino. Capitulo XII	0.5

Samples with higher A-SAS content should be diluted prior to analysis.

The method can be applied to almost any type of water: natural, drinking and wastewater.

METHOD

The fluorimetric method of determination of mass concentration of anionic surfactants in water is based on chloroform extraction of pairs of A-SAS with Acridine Yellow G dye. The intensity of fluorescence of the formed complex is measured by the FLUORAT-02 analyzer and displayed as anionic surfactants concentration in mg/L. The result appears on the PC-operated FLUORATE software.

Samples should be analyzed within 24 hours after taking. Samples prior to analysis should be kept at 4–6°C.

HIGHLIGHTS OF THE FLUORIMETRIC METHOD

- Affordable price for instrument and reagents
- Detection limit is lower and measurement range is wider compared with conventional photometric techniques

EQUIPMENT AND REAGENTS

The following equipment and reagents are used for measuring:

- FLUORAT-02 analyzer with FLUORATE software
- Lumex Instruments optical filters*
- RM of anionic surfactants ion solution (1 g/L)*
- Acridine Yellow G (CAS No. 135-49-9), ≥90%*
- Reagent water complying with grade 1 requirements as defined in ISO 3696
- Chloroform, puriss.
- Hydrochloric acid, puriss.
- Sodium hydroxide, p.a.
- Hexane, puriss.

* - included in Lumex Instruments "Anionic surfactants in water" set, order code 300002513

The contents of this paper are subject to change without notice. To get more specific information, please contact the representative by <u>sales@lumexinstruments.com</u>

IBAEN01.05.03-